

Fast Red TR Salt Tablets (5mg / Tablet)

Catalog No.:	DMR 0440T
Product:	Fast Red TR Salt Tablets
Introduction:	Fast Red Tablets are used as chromogen for alkaline phosphatase (ALP) based immunostaining systems. ALP hydrolyses naphthol phosphate esters to phenolic compounds and phosphates. The phenols couple to colorless diazonium salts to produce insoluble colored azo dye. Fast Red produces a red-colored end product.
Format/Appearance:	20 concave white to light beige 5mg tablets
Composition:	Fast Red TR Salt 5mg Sodium Chloride Remainder
Storage:	Store frozen at -20 to 0°C. Keep dry (suggestion: use dessicant). Product is light sensitive; protect from exposure to light and store in opaque bottle or in dark environment. Warm to room temperature before using.
Precautions and Safety:	Poisonous and suspected carcinogen. Irritant. Avoid contact with skin. Use chemical fume hood with adequate ventilation. Keep away from fire; emits toxic fumes under fire conditions. Follow environmental regulations for disposal. Please contact DBS for MSDS.
Regulatory Information:	
European Information	Caution: Substance not fully tested. Irritant; irritating to eyes, respiratory system, and skin. R: 36/37/38 S: 36/37/39 Wear suitable personal protective equipment, including gloves and eye/face protection.
US Information	This product is subject to SARA Section 313 reporting requirements.

IVD: For In Vitro Diagnostic Use

DBS will not be held responsible for patent infringement or other violation that may occur with the use of our product

DBS

1020 Serpentine Lane, # 114, Pleasanton, CA 94566 Tel: 925 484 3350, Fax: 925 484 3390

Website: www.dbiosys.com e-mail: customersupport@dbiosys.com